

Product Information

MISSION® Lentiviral Packaging Mix

Catalog Number **SHP001**
 Storage Temperature -20 °C

TECHNICAL BULLETIN

Product Description

The MISSION Lentiviral Packaging Mix is an optimized formulation of two plasmids expressing the key HIV packaging genes and a heterologous viral envelope gene.

MISSION lentiviral particles are generated from three main components:

1. The packaging vector, which contains the minimal set of lentiviral genes required to generate the virion structural proteins and packaging functions
2. The vesicular stomatitis virus G-protein (pCMV-VSV-G) envelope vector, which provides the heterologous envelope for pseudo-typing
3. The shRNA transfer vector, which contains the sequence of interest as well as the *cis* acting sequences necessary for RNA production and packaging.

The Lentiviral Packaging Mix contains the first two components; it is designed to be co-transfected along with a compatible lentiviral transfer vector in order to create high-titer pseudo-typed lentiviral particles used for downstream transduction applications.

The MISSION shRNA product line is centered on a transfer vector-based RNAi library against annotated mouse and human genes that can be used in conjunction with this Lentiviral Packaging Mix. In this packaging system, the pseudo-typing with VSV-G broadens the viral tropism associated with the virus.¹ Therefore, these lentiviral particles can efficiently deliver the transfer sequence of interest to a wide variety of cell types, including primary and non-dividing cells.^{2,3}

The particles are produced using a third-generation packaging system, which has many features that lead to enhanced biosafety.⁴ Some of these features are described in Table 1.

Table 1.
Biosafety Features of the Third-Generation Lentiviral Packaging System

Feature	Result
Multi-plasmid approach	No single plasmid contains all the genes necessary to produce packaged lentivirus. Resultant particles are replication-incompetent
Deletion in U3 portion of 3' LTR which eliminates the promoter-enhancer region	Avoids promoter interference issues and further negates the possibility of viral replication
Elimination of the majority of lentiviral genes (Δ vpr, vif, vpu and nef) ⁵	Removes virulence genes which are not necessary for shRNA packaging systems

The MISSION Lentiviral Packaging Mix and the lentiviral shRNA transfer vectors can be easily co-transfected into a mammalian production cell line. We recommend HEK293T cells, ATCC® Number CRL-11268, as they are readily transfected and have been demonstrated to consistently produce a high titer of virus.⁶

Recombinant lentiviruses produced with the MISSION Lentiviral Packaging Mix have not been shown to produce replication competent viral particles because of designed safety features. Users should consult and observe their own institutional guidelines for working with such viral systems.

Reagents

Lentiviral Packaging Mix, 0.25 or 1.7 ml
Catalog Number SHP001

Reagents and equipment required, but not provided

- Lentiviral-based transfer vector(s).
- FuGENE®6 Transfection Reagent, Roche Diagnostics
- HEK293T producer cells
ATCC Number CRL-11268
- Serum-free DME medium
- DME medium supplemented with 10% fetal bovine serum and 4 mM L-glutamine.
- HIV-1 p24 Antigen ELISA Kit, ZeptoMetrix, ZMC Catalog # 0801111

Follow distributor's instructions for culturing and maintaining the producer cell line. It is recommended to use the lowest passage # of cells as possible for transfection experiments.

Precautions and Disclaimer

This product is for R&D use only, not for drug, household, or other uses. Though the lentiviral transduction particles produced are replication incompetent, it is highly recommended that they be treated as **Risk Group Level 2 (RGL-2)** organisms.⁷ Follow all published RGL-2 guidelines for handling and waste decontamination. Also, use extra caution when packaging lentiviral particles that express shRNA-targeting genes involved in cell cycle control, e.g., tumor suppressor genes.

Storage/Stability

If stored properly, shelf life is 2 years from the date of shipment. The Lentiviral Packaging Mix should be stored at -20°C . Avoid multiple freeze thaw cycles.

Procedure Overview

Day 1	Day 2	Day 3	Day 4	Day 5
Plate HEK293T cells	Transfect cells with packaging mix and transfer vector	Re-feed cells with fresh media	First lentiviral particle harvest	Second lentiviral particle harvest; p24 assay

Procedure

Day 1:

Plate the HEK293T cells in complete DME media, supplemented with 10% fetal bovine serum and 4 mM L-glutamine, 24 hours prior to transfection. Cells should reach 50-70% confluency the day of transfection. Recommended cell seeding densities for different plate formats are outlined in Table 2.

Table 2.
Recommended Cell Densities for Various Formats

Format	Initial number of Cells Plated	Final Volume of viral harvest (Assumes two harvests)
96-well plate	16,000 – 20,000 cells/well	200 μL
60 mm dish	5×10^5 cells/dish	8.0 mL
100 mm dish	1×10^6 cells/dish	22 mL
6-well plate	2.3×10^5 cells/well	6.0 mL
T75 flask	1.8×10^6 cells/flask	30 mL
T225 flask	5.4×10^6 cells/flask	90 mL

Day 2:

1. Thaw the vial of Lentiviral Packaging Mix at room temperature. Once thawed, place on ice.

Preparing Volume (in μL) of Lentiviral Packaging Mix

To calculate the amount of Lentiviral Packaging Mix needed for the experiment, refer to Table 3. Prepare excess volume of Lentiviral Packaging Mix based on format used, in order to compensate for pipetting error.

2. Within a sterile polypropylene tube(s) add the calculated volume of Lentiviral Packaging Mix for each transfection.
3. Keep tube(s) on ice.

Table 3.
Volume of Lentiviral Packaging Mix Required for Various Formats

Format	Lentiviral Packaging Mix
96-well plate	1 μL /well
60 mm dish	10 μL /dish
100 mm dish	26 μL /dish
6-well plate	4.6 μL /well
T75 flask	34 μL /flask
T225 flask	107 μL /flask

Preparing Amount (in μg) of Transfer Vector

To calculate the amount of transfer vector needed for the experiment, refer to Table 4. Prepare excess amount of transfer vector based on format used, in order to compensate for pipetting error.

4. Add the calculated volume of transfer vector for each transfection to the sterile polypropylene tube(s).
5. Keep tube(s) on ice.

Table 4.
Amount of Transfer Vector Required for Various Formats

Format	Transfer Vector
96-well plate	0.1 $\mu\text{g}/\text{well}$
60 mm dish	1.0 $\mu\text{g}/\text{dish}$
100 mm dish	2.6 $\mu\text{g}/\text{dish}$
6-well plate	0.5 $\mu\text{g}/\text{well}$
T75 flask	3.4 $\mu\text{g}/\text{flask}$
T225 flask	10.7 $\mu\text{g}/\text{flask}$

Preparing Volume of FuGENE 6 Transfection Reagent/Serum-Free Media Solution

Note: FuGENE 6 transfection reagent has to be prepared in combination with serum-free DME medium

To calculate the amount of FuGENE 6 transfection reagent/serum-free medium, refer to Table 5. Prepare an excess volume based on format used, in order to compensate for pipetting error.

Table 5
Volume of DME Serum-free Medium and FuGENE 6 Transfection Reagent Required for Various Formats

Format	DME Serum-free Medium	FuGENE 6 Transfection Reagent
96-well plate	15 $\mu\text{L}/\text{well}$	0.6 $\mu\text{L}/\text{well}$
60 mm dish	70 $\mu\text{L}/\text{dish}$	6 $\mu\text{L}/\text{dish}$
100 mm dish	182 $\mu\text{L}/\text{dish}$	16 $\mu\text{L}/\text{dish}$
6-well plate	30.3 $\mu\text{L}/\text{well}$	2.7 $\mu\text{L}/\text{well}$
T75 flask	237 $\mu\text{L}/\text{flask}$	21 $\mu\text{L}/\text{flask}$
T225 flask	710 $\mu\text{L}/\text{flask}$	62 $\mu\text{L}/\text{flask}$

6. In a sterile polypropylene tube(s), first add the calculated amount of DME Serum-Free media.

7. To the media, add the FuGENE 6 transfection reagent.
Note: Use polypropylene material tube(s) rather than polystyrene because FuGENE 6 transfection reagent has a tendency to stick to polystyrene.
8. Mix components gently.

Making Transfection Cocktail

Combine all transfection cocktail components together in one polypropylene vial.

9. Add volume of Lentiviral Packaging Mix per one sample based on the recommendation in Table 3.
10. Add amount of transfer vector per one sample based on the recommendation in Table 4
11. Add volume of DME serum-free media/ FuGENE 6 per one sample based on the recommendation in Table 5.
12. Mix gently by pipetting up and down.
13. Incubate transfection cocktail for 15 minutes at room temperature.
14. Add entire transfection cocktail volume from each polypropylene well(s), plate(s), dish(es) and flask(s) to the corresponding well(s), plate(s), dish(es) and flask(s) containing HEK293T cells.

Day 3

Re-feed with fresh media

15. Pre-warm a sufficient amount of complete media to 37 °C to be used for feeding of the cells.
16. At 16 hours post-transfection, remove media from the transfected cells (avoid disturbing cells) and replace with an equal volume of the pre-warmed complete media.
17. Add the volume of DME complete media to each well, dish or flask based on the recommendations in Table 6.

Table 6
Volume of DME Complete Medium Required for Various Formats

Format	DME Complete Medium
96-well plate	100 $\mu\text{L}/\text{well}$
60 mm dish	5 mL/dish
100 mm dish	10 mL/dish
6-well plate	2-3 mL/well
T75 flask	15 mL/flask
T225 flask	45 mL/flask

18. Incubate cells in incubator (37 °C and 5% CO₂) for an additional 24 hours.

Day 4**First harvest of Viral Particles**

19. Pre-warm a sufficient amount of complete media to be used for feeding of the cells.
20. Between 36-48 hours post-transfection, collect viral particles by carefully removing the media and placing it in a collection tube.

The first harvest of viral particles can be stored at 2-8 °C for 24-48 hours. For long-term storage, freeze at -70 °C.

Note: Multiple freeze thaw cycles of viral particles may reduce the infectious viral titer by 20-50% per cycle.

Note: Harvest time points should be optimized depending upon cell health, passage number, and media composition at the time of transfection.

21. Add an equal volume of fresh pre-warmed complete media to cells and incubate for an additional 24 hours

Day 5**Second harvest of Viral Particles**

22. Between 60-72 hours post-transfection collect viral particles by carefully removing the media (if desired, pool with first harvest).

The second harvest of viral particles can be stored at 2-8 °C for 24-48 hours. For long-term storage, freeze at -70 °C.

Note: Multiple freeze thaw cycles of viral particles may reduce the infectious viral titer by 20-50% per cycle.

23. Titer viral particles by performing the HIV p24 Antigen ELISA assay immediately after pooling (see manufacturer's protocol).

References

1. Burns, J.C., *et al.*, Vesicular Stomatitis Virus G Glycoprotein Pseudotyped Retroviral Vectors: Concentration to a Very High Titer and Efficient Gene Transfer into Mammalian and Non-mammalian Cells. *Proc. Natl. Acad. Sci. USA*, **90**, 8033-8037 (1993).
2. Abbas-Terki, T., *et al.*, Lentiviral-mediated RNA interference. *Hum. Gene Ther.*, **13**, 2197-2201 (2002).
3. Stewart, S.A., *et al.*, Lentivirus-delivered stable gene silencing by RNAi in primary cells. *RNA*, **9**, 493-501 (2003).
4. Dull, T., *et al.*, A third generation lentivirus vector with a conditional packaging system. *J. Virol.*, **72**, 8463-8471 (1998).
5. Zufferey, R., *et al.*, *Nat. Biotechnol.*, **15**, 871-875 (1997).
6. Pear, W.S., *et al.*, Production of high-titer helper-free retroviruses by transient transfection. *Proc. Natl. Acad. Sci. USA*, **90**, 8392-8396 (1993).
7. NIH Guidelines for Research Involving Recombinant DNA Molecules (NIH Guidelines) 2002 (<http://www4.od.nih.gov/oba>).
8. Whither RNAi? *Nature Cell Biology*, **5**, 489-490 (2003).

Troubleshooting Guide

Problem	Cause	Solution
Low viral titer or no viral titer determined in HIV p24 Antigen ELISA assay	Lentiviral Packaging Mix was not added to the transfection mix.	Repeat experiment.
	Lentiviral Packaging Mix was not stored properly.	Check if Lentiviral Packaging Mix was stored at –20 °C upon receiving.
	Lentiviral Packaging Mix went through multiple freeze-thaw cycles.	After receiving, generate aliquots of the Lentiviral Packaging Mix and store at –20 °C until ready for use.
	The target gene is essential for cell growth and viability.	Be sure that target gene is not essential for cell growth and viability.
	shRNA construct is larger than 2 kb.	shRNA clones larger than 2 kb can negatively impact production of lentiviral particles.
	Low yield of shRNA plasmid DNA.	The viral-based vectors are known to be difficult to purify. We recommend streaking the bacterial stock on LB/carbenicillin plate to isolate a single colony and DNA purification with GenElute™ HP Midiprep Kit, Catalog Number NA0200, or GenElute HP Maxiprep Kit, Catalog Number NA0300.
	Used poor quality plasmid DNA.	Do not use plasmid DNA from a mini-prep for transfection. Use DNA purification with Sigma's GenElute plasmid purification kits mentioned above
	Seeding density of cells was not optimal.	Follow cell density recommendations in Table 1.
	Media and reagents were stored improperly.	Store media and reagents at temperatures designated by the manufacturers.

Control Selection Table

Sigma's recommended controls for any shRNA experiment are closely aligned with the controls suggested in the *Nature Cell Biology* editorial.⁸

Recommended Control	Objective
Negative Control: Untreated Cells	Untreated cells will monitor cells culture growth conditions and provide a reference point for comparing all other samples.
Negative Control: Transfection with empty pLKO.1 vector, containing no shRNA insert	MISSION pLKO.1-puro Control Vector, Catalog Number SHC001. This vector does not contain an shRNA insert. Particles generated with this construct serve as a useful negative control during transduction. This controls for transduction effects without subsequent activation of the RNAi pathway.
Negative Control: Transduction with Non-targeting shRNA	MISSION Non-Target shRNA Vector Control, Catalog Number SHC002. The vector contains an shRNA insert that does not target any human or mouse genes due to 5 base pair mismatches with known genes in those genomes. This controls for non-specific events during transduction and activation of the RNAi pathway.
Positive Control: Transfection with positive reporter vector	MISSION TurboGFP Control Vector, Catalog Number SHC003. This is a useful positive control for measuring transfection efficiency and optimizing transduction. The TurboGFP Control vector consists of the lentiviral backbone vector, pLKO.1-puro, containing a gene encoding TurboGFP, driven by the CMV promoter. Co-transfection of this control along with the Lentiviral Packaging Mix provides fast visual confirmation of successful transfection and transduction.
Positive Control: Transfection with shRNA targeting reporter vector	MISSION TurboGFP shRNA Control Vector, Catalog Number SHC004. The TurboGFP shRNA vector consists of the pLKO.1-Puro vector, containing shRNA that targets TurboGFP, and can be used as a positive control to quickly visualize knockdown. This TurboGFP shRNA Control Vector has been experimentally shown to reduce GFP expression by 99.6% in HEK 293T cells after 24 hours. Because this vector targets TurboGFP, and it does not target any human or mouse genes, it can also be used as a negative non-target control in shRNA experiments

ATCC is a registered trademark of the American Type Culture Collection.

FuGENE is a registered trademark of Roche Diagnostics

GenElute is a trademark of Sigma-Aldrich Biotechnology

MISSION is a registered trademark of Sigma-Aldrich Biotechnology LP and Sigma-Aldrich Co.

TurboGFP is a trademark of Evrogen Co.

Limited Use Licenses

Sigma has acquired necessary key licenses for lentiviral systems and RNAi and provides freedom to operate under our label license for relevant purchased products. Because Sigma actively evaluates this rapidly evolving intellectual property space, please visit www.sigma.com/shrna for up-to-date information on current licenses for the MISSION® shRNA collections.

Use of this product for Commercial Purposes requires a license from Sigma-Aldrich Corporation. The purchase of this product conveys to the buyer the nontransferable right to use the purchased amount of the product and components of the product in research conducted by the buyer (whether the buyer is an academic or for-profit entity). The buyer cannot sell or otherwise transfer (a) this product (b) its components or (c) materials made using this product or its components to a third party, or otherwise use this product or its components or materials made using this product or its components for Commercial Purposes. Commercial Purposes means any activity by a party for consideration, but excludes not-for-profit core facilities providing services within their own research institutions at cost. Core facilities are invited to join Sigma-Aldrich's RNAi Partnership Program. Details of Sigma-Aldrich's RNAi Partnership Program can be found at www.sigma.com/rpp.

This product is licensed under U.S. Pat. Nos. 5,817,491; 5,591,624; 5,716,832; 6,312,682; 6,669,936; 6,235,522; 6,924,123 and foreign equivalents from Oxford BioMedica (UK) Ltd., Oxford, UK, and is provided for use in academic and commercial *in vitro* and *in vivo* research for elucidating gene function, and for validating potential gene products and pathways for drug discovery and development, but excludes any use of LentiVector® technology for: creating transgenic birds for the purpose of producing useful or valuable proteins in the eggs of such transgenic birds, the delivery of gene therapies, and for commercial production of therapeutic, diagnostic or other commercial products not intended for research use where such products do not consist of or incorporate a lentiviral vector. Information about licenses for commercial uses excluded under this license is available from Oxford BioMedica (UK), Ltd., Medawar Center, Oxford Science Park, Oxford OX4 4GA UK enquiries@oxfordbiomedica.co.uk or BioMedica Inc., 11622 El Camino Real #100, San Diego CA 92130-2049 USA. LentiVector is a registered US and European Community trademark of Oxford BioMedica plc.

This product (based upon the lentikat system) is sub-licensed from Invitrogen Corporation under U.S. Patent Nos. 5,686,279, 5,834,256, 5,858,740; 5,994,136; 6,013,516; 6,051,427, 6,165,782, and 6,218,187 and corresponding patents and applications in other countries for internal research purposes only. Use of this technology for gene therapy applications or bioprocessing other than for nonhuman research use requires a license from Cell Genesys, Inc. Please contact Cell Genesys, Inc. at 342 Lakeside Drive, Foster City, California 94404. Use of this technology to make or sell products or offer services for consideration in the research market requires a license from Invitrogen Corporation, 1600 Faraday Ave., Carlsbad, CA 92008.

RC,SG,SU,PHC 11/12-1